


# MixAmatic

by Pepetools

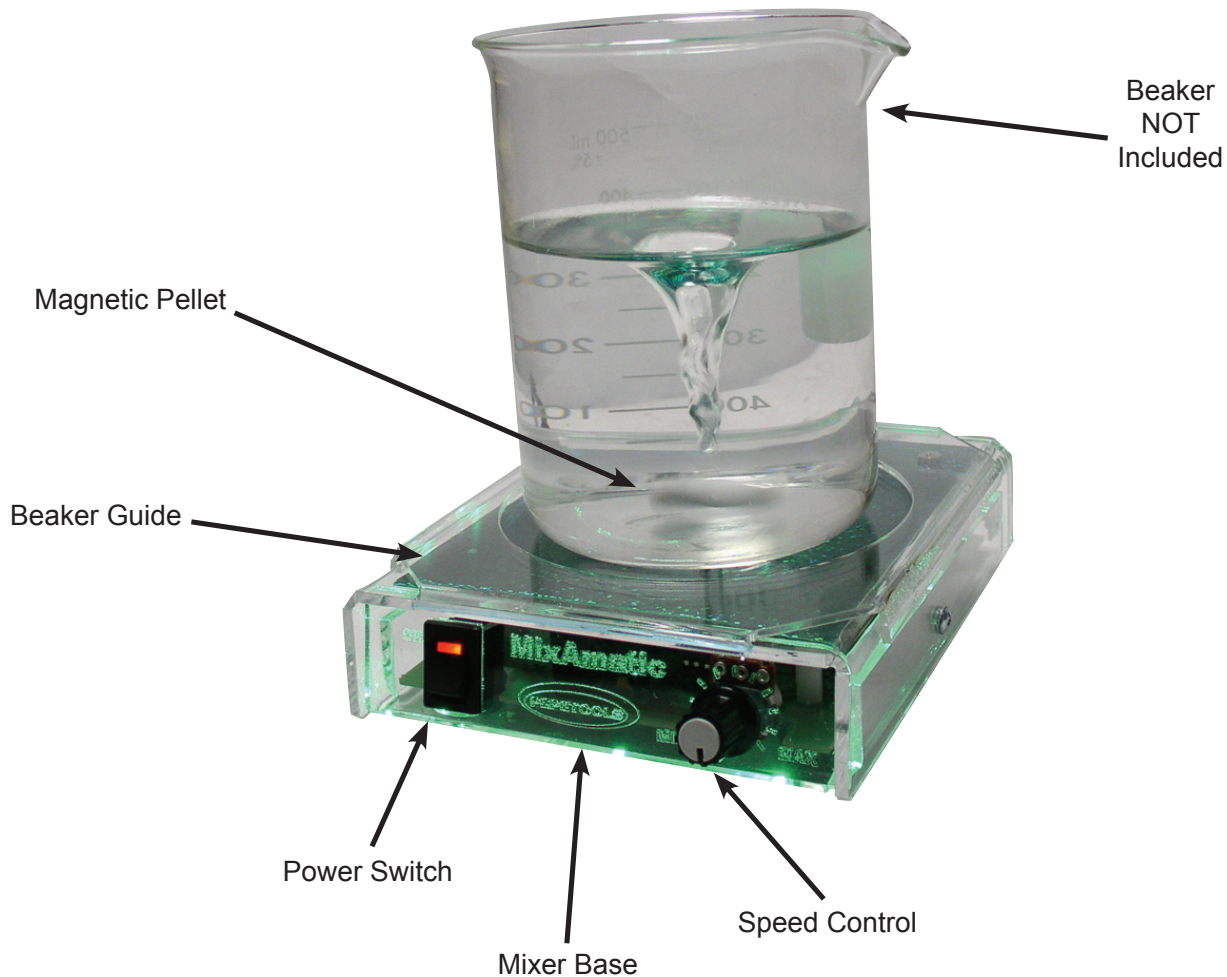

## Setup:

This unit is designed to be used with a 1000ml beaker., but will work well with smaller sizes as well.

Begin by placing the beaker guide and pad on the top of the Mixer. Place the magnetic Pellet into the center of the beaker . Fill your beaker with your solution. Plug in your power cord to the back of the unit.

## Use:

To use the mixer first ensure that the speed control knob is turned all the way to the left. Turn on the power and turn the knob to the right to increase the speed to your desired level of agitation. Ideally, the vortex created by the mixer should not be very deep, a vortex that barely breaks the liquid's surface works best.

## Tip:

If the pellet moves out of the center of the beaker, simply turn the speed all the way down, then increase it slowly and the pellet will be pulled back to center.